



T-Blocks

Product information

Tissue Blocks

Tissue Blocks (T-Blocks) are porous hydrogel scaffolds with polarized microstructures that enable primary cells to attach, migrate, and proliferate after seeding. As primary cells require more surface area to grow, additional T-Blocks may be joined to the original T-Block seeded with cells. Cells will migrate and expand within additionally joined T-Blocks. T-Blocks can be used to maintain primary cell cultures without the need to sub-culture when cells become confluent. See **Table 1** for instructions on how to properly store T-Blocks.

Table 1. Contents and storage

Material	Amount	Storage	Stability
PEGDA	1 vial containing 1 T-Block (1 g)	<ul style="list-style-type: none">• 9 – 30 °C, Max 5-day storage• 4 - 8 °C, long-term storage• Sterile Distilled H₂O	When stored as directed, product is stable for 6 months
* Note that vial contents are hardly visible.			

Properties of T-Blocks

T-Blocks are porous hydrogel scaffolds that enable efficient growth of primary cells in 3D. T-Blocks are made out of biocompatible hydrogel compositions which include proprietary blends of polyethylene glycol diacrylate (PEGDA). T-Blocks coated with human extracellular matrix coatings facilitate attachment and migration of primary cell populations. T-Blocks are a platform by which the T-Block substrate can be expanded by joining additional T-Blocks to the original T-Block. This process facilitates migration and proliferation of primary cells, while maintaining the primary cell phenotype and eliminating the requirement of sub-culture.

Materials required, but not provided.

- Sterile Hanks Balanced Salt Solution (HBSS) [without Ca²⁺ and without Mg²⁺], pH 7.2
- Human Fibronectin (Corning Cat. No. 354008)
- Primary cell population of choice
- Sterile Expansion media of choice
 - Note: Use serum-free media only, if cells are intended to be removed from T-Block
- Sterile Trypsin, 0.05% EDTA
- Sterile spatula
- Single channel pipettors



For Research Use Only. Not for human or diagnostic use.
Publication Number: T-0001 Ver. A.011

- Pipettor tips
- 50-mL Conical Tubes

Methods

Perform all procedures using aseptic technique unless otherwise noted.

Pre-Soak T-Blocks

- Submerge T-Blocks in Hanks Balanced Salt Solution (HBSS) [without Ca^{2+} and without Mg^{2+}], and soak T-Blocks one day prior to coating.



Coating T-Blocks with Fibronectin

- Equalize Fibronectin to room temperature before use.
- Reconstitute fibronectin solution to a stock concentration of 1 mg/mL in HBSS [without Ca^{2+} and without Mg^{2+}].
 - o Gently swirl the solution for 1 minute to mix.
 - DO NOT VORTEX
- Create a 50 $\mu\text{g/mL}$ working fibronectin solution in HBSS [without Ca^{2+} and without Mg^{2+}].
 - o Gently swirl the solution for 1 minute.
- Place T-Block in working fibronectin solution for at least 6 hours.
 - o Make sure the T-Block is completely submerged.
 - o T-Blocks may be incubated with fibronectin overnight.
 - o T-Blocks may be coated in polypropylene conical tubes instead of well plates or mono plates.
- Remove T-Block from working fibronectin solution with sterile spatula.
- Gently wash the T-Block with sterile HBSS [without Ca^{2+} and without Mg^{2+}] three times, before seeding cells.


Seeding T-Blocks with Primary Cells

- Place coated T-Blocks in sterile non-tissue culture treated 6-well plate with the T-Block bottom “rough” side facing plate floor.
- Suspend primary cells of choice in media of choice at a concentration of 100,000 cells to 1,000,000 cells per 0.5 mL, dependent on cell growth rate.
 - o Use serum-free media exclusively, if cells will be removed from T-Block.
- Gently pipette cell solution onto TOP side of T-Block dropwise.
 - o Top of T-Block has a smooth surface (See **Table 2A**).
 - o Bottom of T-Block has a “rough” appearance. (See **Table 2B**).

Table 2. Top and Bottom of T-Block after Submerged in Media

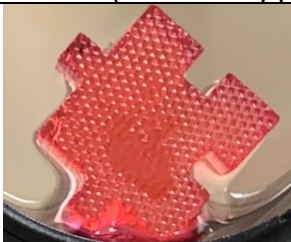A. Top of T-Block (Smooth Appearance)	B. Bottom of T-Block (Rough Appearance)
	

- Wait 5 minutes.
- Add 6.5 - 8.5 mL of media of choice to well containing T-Block.
 - Add media against wall of well. **DO NOT ADD** media to top of T-Block after cells have been added.
 - Ensure that T-Block is fully submerged.
- Check T-Blocks for confluency using an inverted microscope and change media every 2 days.
 - When changing media take care not to disturb the T-Block. For best results, remove media manually with a pipette.
 - If using an aspirating system that utilizes a vacuum pump, be sure to aspirate media as far away from the T-Block as possible. **DO NOT** let T-Block touch aspirating tip or T-Block may be damaged.
- When primary cells occupy approximately 60% of all cell surfaces within T-Block, join an additional T-Block to the original T-Block by sliding the tongue of new T-Block into a groove of original T-Block as shown in **Figure 1**.
 - If needed, carefully transfer T-Block to mono-plate before adding new T-Block to provide more cell culture volume.
 - **ALWAYS** ensure at least one side of each T-Block is exposed to media when joining T-Blocks together to permit proper media diffusion.

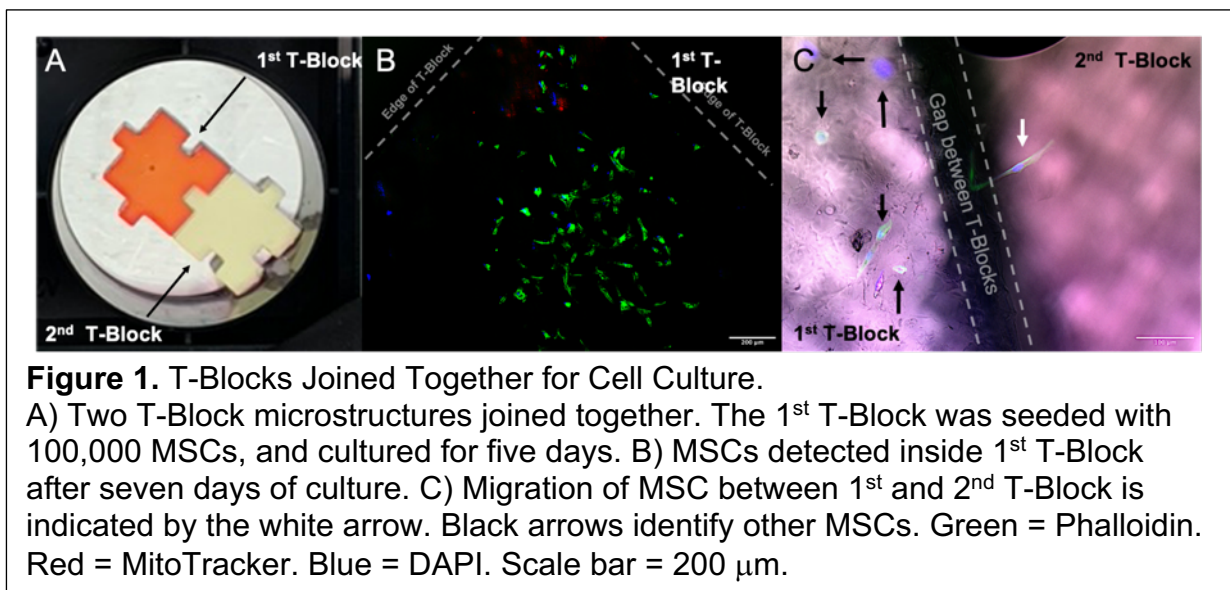

- Continue culturing T-Blocks with preferred media.
 - Ensure that enough media is added to fully submerge T-Block(s).
- Repeat process until desired cell concentration is reached.
- Follow standard trypsinization protocols for dissociating primary cells from T-Blocks when cells are needed for analysis.


- Dispose of used T-Block following standard biological hazardous wastes protocols mandated by institution, locality, state, and federal guidelines.

Supplemental Information

Cell Growth and Attachment

When T-Blocks are properly coated with extracellular matrix and seeded properly, cells will infiltrate pores and migrate throughout the T-Block. Use 10X objectives or greater when viewing cells with bright-field or epi-fluorescent microscopes to minimize light reflection within the T-Block. Cells can be identified by their “rough” texture. Use of fluorescent dyes such as Hoechst 33342 will enhance visualization of cells within T-Block dramatically. **Table 3** illustrates both transmission light with fluorescent overlay and fluorescence markers (Green = Phalloidin, Red = MitoTracker, Blue = DAPI) without transmission light. When viewing cells, keep in mind that all boundaries of cell may not be completely in focus if cell is residing in multiple Z-planes. Adjustment of the fine focus knob on the microscope may help in viewing entire cell body within T-Block.

Table 3. Visualization of T-Blocks with Cells

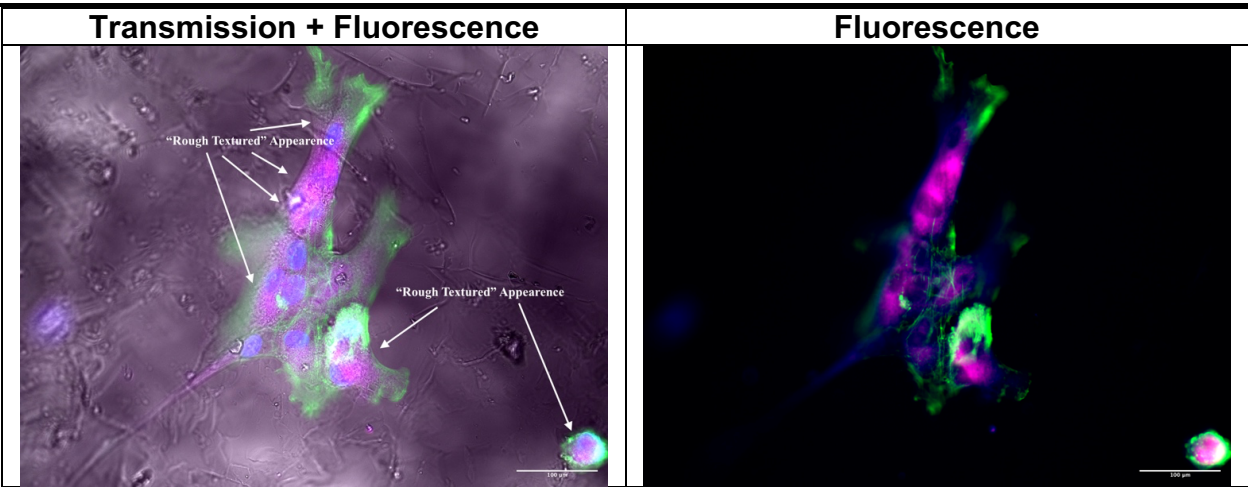



Ordering Information

Cat. No.	Product Name	Unit Size
PAT001	PEG T-Block.....	1 U
PAT002	PEG T-Block.....	4 U
PAT003	PEG T-Block.....	12 U
PAT004	PEG T-Block.....	24 U

Documentation and Support

Customer and Technical Support

Visit ronawk.com for the latest in services and support, including:

Contact telephone numbers

Product support, including:

- Product FAQs
- Training

Order and web support

Product documentation, including:

- User guides, manuals, and protocols
- Certificate of Analysis
- Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited Product Warranty

Ronawk LLC and/or its affiliate(s) warrant their product as set forth in the Ronawk LLC General Terms and Conditions of Sale found on the Ronawk LLC website at www.ronawk.com/products/termsandconditions.html. If you have any questions, please contact Ronawk LLC at ronawk.com/product/t-block/.

The information in this guide is subject to change without notice.

For Research Use Only. Not for human or diagnostic use.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, RONAWK LLC AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Manufacturer: Ronawk LLC | 500 N. Rogers Road, Suite 500-B20 | Olathe, KS 66062

Trademarks: All trademarks are the property of Ronawk LLC and its subsidiaries unless otherwise specified.

©2020 Ronawk LLC. All rights reserved.